




**Thermo Scientific
Environmental Chambers**
Controlled Temperature Storage Solutions

Optimum uniformity
plus a wide temperature range

Thermo
SCIENTIFIC

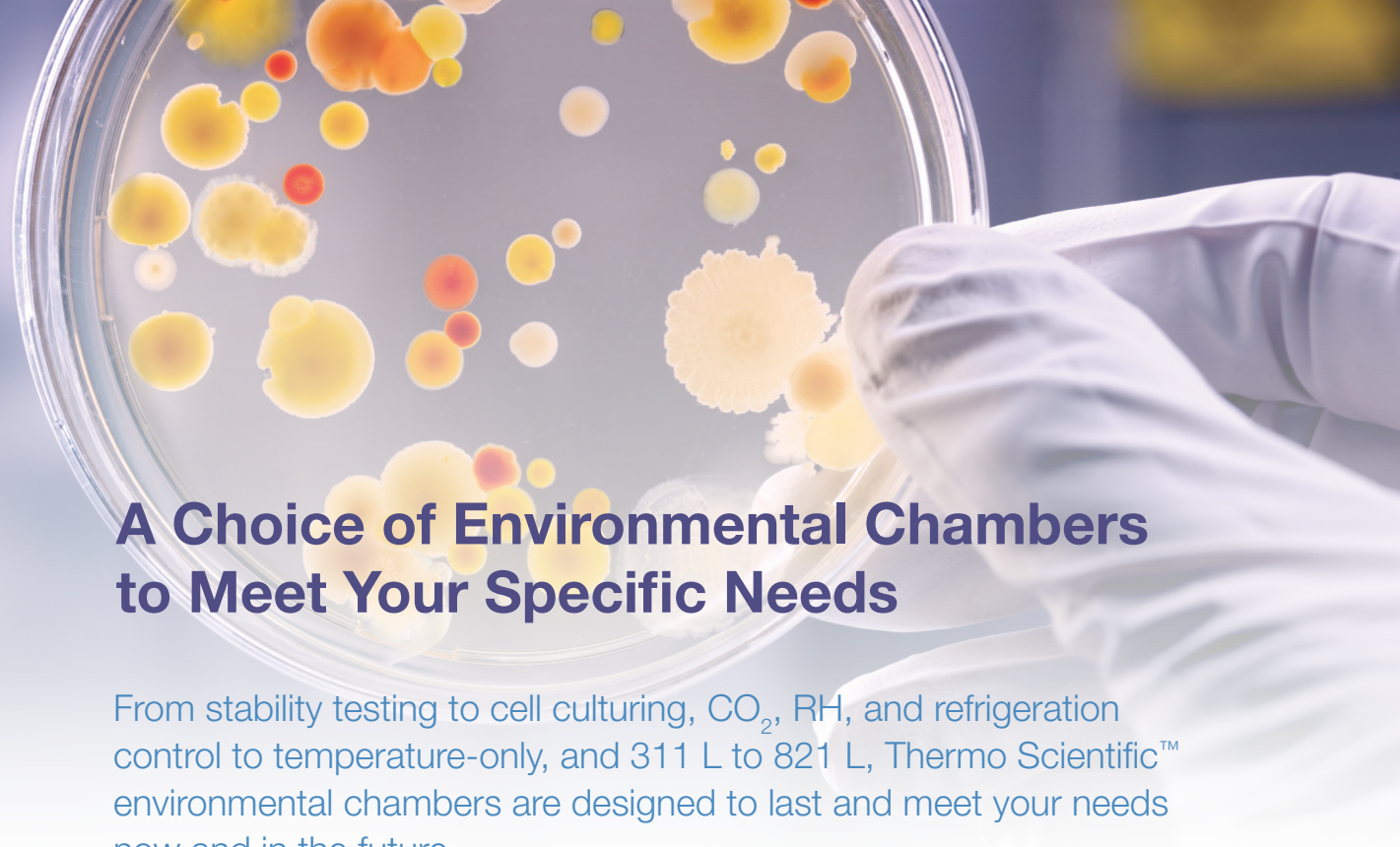

A Choice of Environmental Chambers to Meet Your Specific Needs

From stability testing to cell culturing, CO₂, RH, and refrigeration control to temperature-only, and 311 L to 821 L, Thermo Scientific™ environmental chambers are designed to last and meet your needs now and in the future.

If you require a high precision cabinet to meet ICH, FDA, TAPPI, DIN12880 or national testing standards, our **Thermo Scientific 3948/3907 environmental chamber series** is the ideal choice.

- Optimum uniformity and a wide temperature range designed for easy compliance with regulatory requirements.
- Rugged construction and durable shelving for long life.
- Broad range of options to create the perfect fit for your applications.

Consider our **Thermo Scientific 3962 reach-in incubator/environmental chamber** if you are focused on elevated temperature-only segments of drug stability and culture.

- Optimum uniformity and an elevated temperature range designed for easy compliance with regulatory requirements.
- Rugged construction and durable shelving for long life.
- Valuable features for flexibility and cost-effectiveness.

At-A-Glance Chamber Comparison

| Type | Thermo Scientific Cat. No. | Temp Range | Uniformity | RH | Size (L) | CO ₂ |
|-----------|----------------------------|---------------------------|---|-----------------------------|----------|-----------------|
| Stability | 3948 | 0°C TO 60°C | ±0.3°C @ 25°C to 37°C | Above ambient to 95% @ 37°C | 821 | Optional |
| | 3907 | 0°C TO 60°C | ±0.3°C @ 25°C to 37°C | Above ambient to 95% @ 37°C | 311 | Optional |
| | 3962 | 5°C above ambient to 60°C | ±0.4°C @ 30°C ±0.3°C @ 37°C ±0.5°C @ 45°C | N/A | 821 | N/A |

Performance values are based on 15 point measurement at ambient temperature of 22°C and 230V line voltage.

Thermo Scientific Stability Chambers for a Multitude of Applications

This series of environmental chambers meets International Conference on Harmonization (ICH) guidelines for drug substance and product storage testing. The units are ideal for drug stability studies, shelf life testing for packaged products, and insect and other large-scale biological research.

Optimum uniformity and a wide temperature range designed for easy compliance with regulatory requirements

These chambers feature a 0°C to 60°C (32°F to 140°F) temperature range that can be set in 0.1°C increments. The broad temperature range, and optimum temperature uniformity and recovery contribute to an ideal test environment, even when the chamber contains large product loads and low output, heat-generating equipment.

Advantages of our directed horizontal laminar airflow system

Our directed airflow system (shown below) promotes an ideal growth environment. The enhanced design includes a positive pressure feed


plenum on the right side of the chamber and a negative pressure return plenum on the left to distribute the airflow uniformly throughout the chamber.

This combination directs air across the surface of each solid shelf. Even when filled with samples or equipment, each shelf receives a consistent flow of conditioned air for optimum temperature uniformity and recovery.

As opposed to our horizontal airflow system, top-to-bottom (non-directed) airflow systems use a top-mounted fan to push air down through wire shelves.


Temperature uniformity and recovery can deteriorate quickly when shelves are filled because air movement is blocked. That variation in temperature, alone or when combined with frequent door openings, may compromise environmental conditions or make process validation difficult.

Standard 4 to 20 milliamp output connects to most alarm/monitoring systems, allowing you to meet internal and regulatory product documentation requirements. Chart recorders are also available.

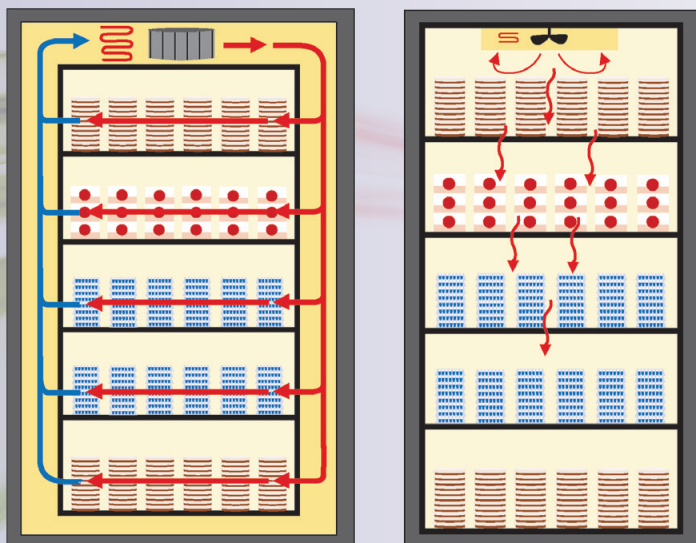

The Thermo Scientific directed airflow design (left) versus a top-to-bottom, undirected airflow design (right).

The directed airflow minimizes the risk of product desiccation and loss, and wasted time and money due to poor temperature uniformity and recovery.

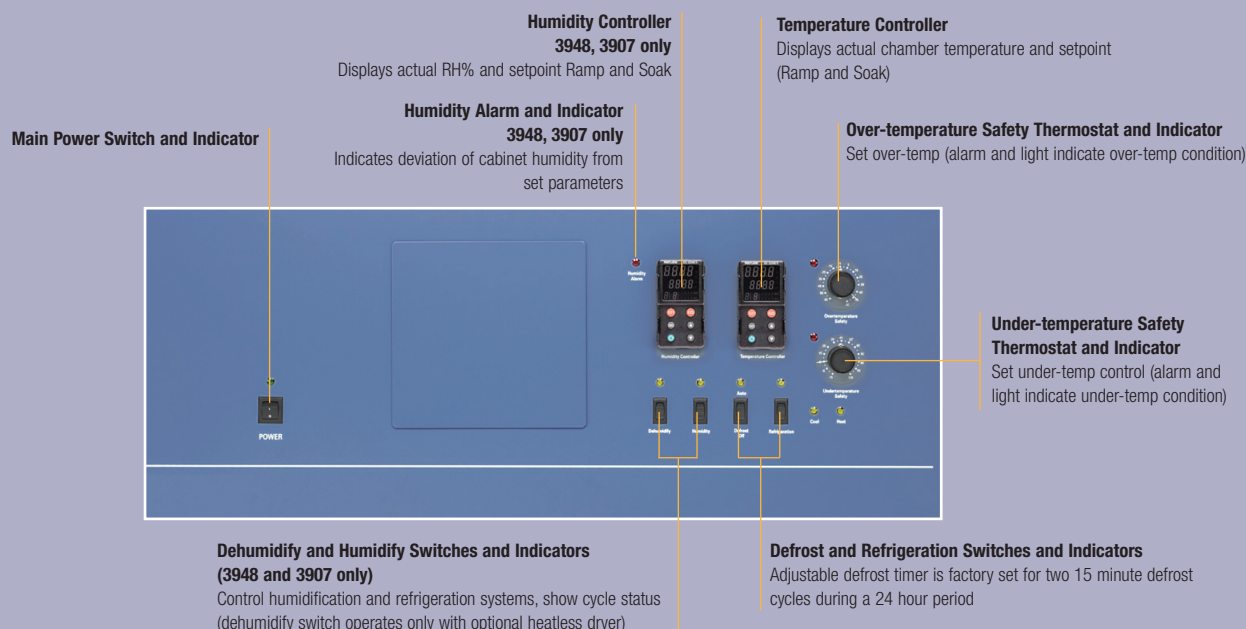

Rugged construction and durable shelving for long life

- Exterior cabinet is 18 gauge cold-rolled steel, powder coated for durability and a high quality appearance that resists scratches and chipping.
- 51 mm (2.0") thick, non-CFC, foamed-in-place polyurethane cabinet insulation provides maximum product thermal protection and increased rigidity for reliable daily use.
- Phenolic-coated evaporator protects against corrosive by-products or agents in the chamber atmosphere (e.g., secretions from insects such as flies), designed for long life.
- Solid, stainless steel interior is more durable and corrosion-resistant than plastic or painted metal, designed for long life and minimizing equipment costs.
- Heavy-duty, solid, stainless steel shelves are easy-to-clean, saving time and effort; more corrosion-resistant than coated wire shelves for long life; and adjustable for convenience.
- 821 L models include 6 shelves that are adjustable on 76mm (3 in.) centers.
- 311 L model includes 3 shelves that are adjustable on 51mm (2 in.) centers.
- Reinforced floor/shelf components for 821 L models are available so you can use shakers, cell rollers, or other heavy equipment or product loads inside the chamber to meet a variety of applications.
- Swivel, locking casters designed for easy mobility during installation and cleaning. Leveling feet provide stability and added safety in the lab.

Broad range of options to create the perfect fit for your applications

Many options are available so you can tailor this series to your specific applications.

Factory-installed options (ordered separately) include:

- Choice of access ports for inserting data logging sensors or cables.
- Convenience receptacles to run electrical devices in chamber.
- Choice of chart recorders for data logging.
- Choice of data outputs for monitoring/alarm (3960/3962 models only).
- Infrared CO₂ Control Package for control of a CO₂ environment.
- Heatless dryer for low humidity conditions – check temperature / humidity chart (see page 5 for details).
- Stainless steel glass door cover for light sensitive applications.
- Door options include door lock, left-hand door, and Lexan™ inner door.
- Choice of shaker support systems provides system safety.
- Temperature mapping and calibration certificate – customized.

Accessories/on-site options (ordered separately) include:

- Extra and special shelving.
- Reinforced floor options for heavy weights – up to 68 kg (150 lb.).
- Sensaphone™ telephone dialing systems – interfaces with standard touch-tone phone.


Options for the humidity system:

- Condensate pump kit – recommended when the accessible drain on lab wall is well above the drain on back of environmental chamber.
- Condensate evaporator – external heating device which burns off water that drips out the back of the unit when there is no drain available.
- Cell roll system for production of monolayer cell cultures (see page 7 for details).
- On site IQ/OQ/PQ validation services.

See page 11
for order numbers,
model capabilities and
additional information.

Temperature/Humidity Performance Chart for 3948 and 3907

This chart illustrates the maximum and minimum attainable humidity within the cabinet at a given set temperature under the following nominal conditions: refrigeration on, defrost off, no wet bulb recorder, no air exchange, set for 22.2°C (72°F) and 50% RH ambient. Optional heatless dryer is available for lower humidity levels.

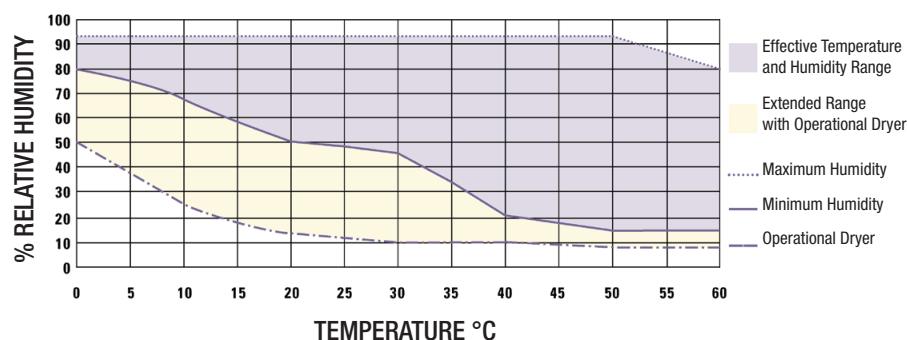



Chamber Specifications

| Thermo Scientific Cat. No. | 3948 ** | 3907** |
|----------------------------|--|--|
| Description | 821 L Incubated/Refrigerated Environmental Chamber with humidity control | 311 L Incubated/Refrigerated Environmental Chamber with humidity control |
| Voltage | 230V, 50/60 Hz | 230V, 50/60 Hz |
| Temperature | | |
| Control | ±0.1°C @ 25°C to 37°C (77°F to 98.6°F) | ±0.1°C @ 25°C to 37°C (77°F to 98.6°F) |
| Range | 0°C to 60°C (32°F to 140°F) | 0°C to 60°C (32°F to 140°F) |
| Sensor | RTD | RTD |
| Controller | Microprocessor, PID | Microprocessor, PID |
| Programmability | Ramp & Soak (to 40 steps) | Ramp & Soak (to 40 steps) |
| Setpoint | Digital | Digital |
| Display | Digital dual LED | Digital dual LED |
| Readability & Setability | 0.1°C | 0.1°C |
| Uniformity* | ±0.3°C @ 25°C to 37°C (77°F to 98.6°F) | ±0.3°C @ 25°C to 37°C (77°F to 98.6°F) |
| Temperature Safety | | |
| Sensor & Controller | Thermostat | Thermostat |
| Setpoint | Analog reference dial | Analog reference dial |
| Alarm | Audible/visual | Audible/visual |
| Refrigeration | | |
| Compressor | 185W (1/4 HP), air-cooled | 185W (1/4 HP), air-cooled |
| Refrigerant | Non-CFC, R134A refrigerant | Non-CFC, R134A refrigerant |
| Humidity | | |
| Input Water Quality | 50K-1Megohm resistance | 50K-1Megohm resistance |
| Input Water Volume | 3.78 L/hour (1 gallon/hour) | 3.78 L/hour (1 gallon/hour) |
| Input Water Pressure | Gravity to 2.76 bar (40 psi) | Gravity to 2.76 bar (40 psi) |
| RH | Above ambient to 95% @ 37°C (98.6°F) ±5.0% RH | Above ambient to 95% @ 37°C (98.6°F) ±5.0% RH |
| Sensor | Bulk polymer | Bulk polymer |
| Controller | Microprocessor, PID | Microprocessor, PID |
| Programmability | Ramp & Soak (to 40 steps) | Ramp & Soak (to 40 steps) |
| Setpoint | Digital (direct set in % RH) | Digital (direct set in % RH) |
| Display | Digital, dual LED (direct read in % RH) | Digital, dual LED (direct read in % RH) |
| Readability & Setability | 1.0% | 1.0% |
| Steam Generator | Initial fill approx. 1 liter (1 qt.) | Initial fill approx. 1 liter (1 qt.) |
| Setable Alarm | Audible/visual | Audible/visual |
| Certification | CE | CE |

*Six shelves in 3948; three shelves in 3907; uniformity is ±0.5°C at all other temperature parameters. Values based on 15 point measurement at ambient temperature of 22°C, 230V line voltage.

**All models require a drain unless fitted with the optional Condensate Evaporator (No. 1900031). They also require a constant water source.

Chamber Specifications continued

| Thermo Scientific Cat. No. | 3948 ** | 3907** |
|------------------------------------|---|---|
| Description | 821 L Incubated/Refrigerated Environmental Chamber with humidity control | 311 L Incubated/Refrigerated Environmental Chamber with humidity control |
| Fittings** | | |
| Fill Port | 1/8" female NPT | 1/8" female NPT |
| Drain Port (bottom) | 3/8" FPT & 3/8" O.D. P-trap right rear | 3/8" FPT & 3/8" O.D. P-trap right rear |
| Unit Heat Load | | |
| 220V | 1750W (6000 BTU/hour) | 1750W (6000 BTU/hour) |
| Dimensions | | |
| Exterior | | |
| Width | 96.5cm (38 in.) | 96.5cm (38 in.) |
| Height | 224.8cm (88.5 in.) | 130.8cm (51.5 in.) |
| F-B | 81.3cm (32.0 in.) | 81.3cm (32.0 in.) |
| Interior | | |
| Width | 78.7cm (31.0 in.) | 78.7cm (31.0 in.) |
| Height | 152.4cm (60.0 in.) | 60.9cm (24.0 in.) |
| F-B | 68.6cm (27.0 in.) | 68.6cm (27.0 in.) |
| Shelves | | |
| Dimensions | 77.8cm x 65.6cm F-B (30.6 in.W x 25.8 in.) | 77.8cm x 65.6cm F-B (30.6 in.W x 25.8 in.) |
| Construction | Solid stainless steel, reinforced | Solid stainless steel, reinforced |
| Surface Area per Shelf | 0.5 sq. m (5.4 sq. ft.) | 0.5 sq. m (5.4 sq. ft.) |
| Max. Surface Area per Chamber | 9.7 sq. m (104.3 sq. ft.) | 5.5 sq. m (59.4 sq. ft.) |
| Standard/Maximum Number of Shelves | 6, 19 | 3, 11 |
| Load per Shelf | 15.9 kg (35 lbs.), slide in/out 22.7 kg (50 lbs.), stationary | 15.9 kg (35 lbs.), slide in/out 22.7 kg (50 lbs.), stationary |
| Construction | | |
| Interior Volume | 821.2 liters / 29.0 cu. ft. | 311.5 liters / 11.0 cu. ft. |
| Interior Material | EN standard 1.4301, 2B, Stainless Steel | EN standard 1.4301, 2B, Stainless Steel |
| Exterior Material | Cold-rolled steel, powder coated | Cold-rolled steel, powder coated |
| Outer Door Gasket | Four-sided, vinyl compression | Four-sided, vinyl compression |
| Electrical | | |
| | 230V,50/60 Hz 15.0 Amps | 230V,50/60 Hz 15.0 Amps |
| Plug | CEE 7/7  | CEE 7/7  |
| Dry Contacts | Common, NO, NC | Common, NO, NC |
| Data Output | 4-20 milliamp, temperature and RH | 4-20 milliamp, temperature and RH |
| Power Switch | 2 Pole | 2 Pole |
| Weight | | |
| Net | 347.0 kg (765 lbs.) | 260.8 kg (575 lbs.) |
| Shipping (motor) | 426.4 kg (940 lbs.) | 324.3 kg (715 lbs.) |

**All models require a drain unless fitted with the optional Condensate Evaporator (No. 1900031) and a constant water source.

Incubator/Environmental Chamber for Elevated Temperature Applications

The 821 L (29 cu. ft.) chamber 3962 offers precise temperature control in a range of 5°C above ambient to 60°C (140°F), making this unit ideal for clinical applications, elevating temperature stability

- Large size for high volumes and a broad range of products.
- Heated glass door for minimal condensation and a clear view.
- Interior and exterior convenience receptacles, and five adjustable shelves.
- Thermo Scientific™ Enviro-Scan™ controls with easy-to-read display of actual temp studies, and shelf life testing.

Model Number 3962



Cell Roll System

The optional cell roll system allows extensive production of monolayer cell cultures in standard roller culture vessels. Oxygenation and exposure of the cells to the media growth area are improved. Culture yields are increased by the uniform temperature control and cell roll system's continuous, gentle rotation.

The environmental chamber accommodates a cell roller up to 7 decks high with 5 positions per deck for a maximum total of 35 positions, or bottles.

Achieving maximum capacity requires option 4862: a three-tier cell roller base (15 positions), four add-on tiers (20 positions), and a reinforced floor/ramp.

All position drive is standard. Adjustable speed control provides precise speeds of 0.125 to 6.25 RPM with $\pm 1.0\%$ accuracy, based on 110mm bottles.

Optimum Uniformity and an Elevated Temperature Range Designed for Easy Compliance with Regulatory Requirements

Our directed airflow system promotes an ideal growth environment. The proven design includes a positive pressure feed plenum and a negative pressure return plenum. This combination directs air across the surface of each solid shelf. Even when filled with samples or equipment, each shelf receives a consistent flow of conditioned air for optimum temperature uniformity and recovery. By design, the feed plenum cannot be blocked by the chamber's contents. (Refer to the diagram on page 3.)

Intuitive, powerful Enviro-Scan



- > Designed to ensure precise temperature control without complicated programming
- > Provides audible/visual under-temperature and over-temperature alarms for peace-of-mind
- > Includes an easy-to-read display for convenient, continuous monitoring
- Enviro-Scan operating modes include Run, Setpoint, Calibration, and System Configuration.
- Standard remote alarm contacts and available data outputs allow connection to an in-house monitor/alarm system to track chamber conditions, helping you meet internal and regulatory documentation requirements. Chart recorders are available.



Durable construction and shelving for long life


- Exterior cabinet is 18 gauge cold-rolled steel, powder coated for durability and a high quality appearance that resists scratches and chipping.
- Maximum product thermal protection and added cabinet strength for reliable daily use are provided by the 51mm (2 in.) thick fiberglass cabinet insulation.
- Solid, stainless steel interior is more durable and corrosion-resistant than plastic or painted metal, ensuring long life and minimizing equipment costs.
- Five heavy-duty, solid, stainless steel shelves, adjustable on 51mm (2 in.) centers, are easy-to-clean, saving time and effort; more corrosion-resistant than coated wire shelves for long life; and adjustable for convenience.
- Reinforced floor/shelf components are available so you can use shakers, cell rollers, or other heavy equipment or product loads inside the chamber to meet a variety of applications.

Valuable features included for cost-effectiveness and flexibility

Casters, access ports, convenience receptacles, heated glass door, and remote alarm contacts – all are included up-front. Our 3962 Incubator/Environmental Chamber is prepared to meet your changing application needs without the added cost of expensive add-on options.

- Swivel, locking casters ensure easy mobility during installation and cleaning. Leveling feet provide stability and added safety in the lab.
 - Thru-wall access ports, located on the right and left sides of the chamber, make it possible to add probes, sensors, etc. without altering the cabinet.
 - Interior and exterior accessory receptacles provide a convenient power source.
 - > Interior receptacle European 230V CEE 7 on 3962 located in the upper right corner of the rear wall, permits the use of shakers, cell rollers, and other equipment inside the chamber, eliminating the inconvenience of an extra external power strip.
 - > Exterior receptacle, located on the upper-right side of the control panel, is available for connecting an optional recorder or other equipment.
- 

Chamber Specifications continued

| | |
|-------------------------------------|---|
| Thermo Scientific Cat. No. | 3962 |
| Description | 821 L Incubated Environmental Chamber |
| Voltage | 200-230V, 50/60 Hz |
| Temperature | |
| Control | ±0.1°C |
| Range | 5°C above ambient to 60°C (140°F) |
| Uniformity | ±0.4°C @ 30°C (86°F) / ±0.3°C @ 37°C (98.6°F) / ±0.5°C @ 45°C (113°F) |
| Low Alarm | User programmable and indicator |
| Over-temperature | |
| Sensor | Precision thermistor |
| Setability | 0.1°C |
| Function | User programmable, action (shuts off heat), and indicator |
| Temperature Safety (Cabinet) | |
| Sensor | Independent thermostat |
| Type | Independent analog electronic |
| Fittings | |
| Access Port With Stopper | 6.1 cm (2.4 in.) I.D., one on each side |
| Unit Heat Load | |
| 115V/230V | 150W (510 BTU / hour); 129 kcal per hour |
| Shelves | |
| Dimensions | 77.7cm x 65.5cm F-B (30.6 in. W x 25.8 in.) |
| Construction | 2B finish, solid stainless steel |
| Surface Area | 0.5 sq. m (5.4 sq. ft.) per shelf |
| Max. per Chamber | 13.0 sq. m (140.4 sq. ft.) |
| Standard/Maximum No. of Shelves | 5, 26 |
| Loading | 13.6 kg (30 lbs.) per shelf, fully inserted and stationary; not to exceed 136.0 kg (300 lbs.) total per cabinet |
| Construction | |
| Interior Volume | 821.2 liters (29 cu. ft.) |
| Interior Material | Type 304, 2B finish, stainless steel |
| Exterior Material | Cold-rolled steel, powder coated |
| Insulation | 5.1cm (2.0 in.) fiberglass |
| Exterior Door | Heated, triple pane, tempered glass |
| Outer Door Gasket | Molded vinyl |
| Electrical | |
| | 200/230V, 50/60 Hz, 5.0 Amps (Operating range 180-250V includes voltage fluctuations) |
| Circuit Breaker/Power Switch | 8 Amps; 2 pole |
| Exterior Convenience Receptacle | 75 Watts maximum (matches cabinet voltage) |
| Interior Convenience Receptacle | 230 Watts maximum (matches cabinet voltage) |
| Plug | Plug 230V: CEE 7/7 Plug  |
| Alarm Contacts | Temperature deviation and power failure; Common, NO, and NC; customer connections through RJ11 jack on rear of unit |
| Data Outputs (opt.) | RS-485, 0-1V, 0-5V, 4-20 milliamp |
| Dimensions | |
| Exterior | 96.5cm x 203.2cm x 83.8cm F-B (38.0 in. W x 80.0 in.H x 33.0 in.) |
| Interior | 78.7cm x 152.4cm x 68.6cm F-B (31.0 in. W x 60.0 in.H x 27.0 in.) |
| Weight | |
| Net | 226.8 kg (500 lbs.) |
| Shipping (motor) | 299.4 kg (660 lbs.) |
| Certification | CE |

Specifications are based on nominal voltages of 230V in ambients of 22°C to 25°C.
Performance is plus or minus the least significant digit unless otherwise specified.

Accessories (All are customer installed unless indicated otherwise)

| Thermo Scientific Cat. No | Description |
|---|---|
| Shelving | |
| Stainless Steel Shelves for 3948, 3907, and 3962 | |
| 224139 | Solid Stainless Steel Shelf and Channels for 3948, 3907 and 3962 |
| 224155 | Perforated Stainless Steel Shelf and Channels for 3948, 3907 and 3962 |
| 224161 | Reinforced Stainless Steel Shelf and Channels for 3948, 3907, and 3962; two per unit max., not for shakers; 68.0 kg load maximum with shelf fully inserted and stationary |
| Shaker Support System, factory installed | |
| 1900005* | Two shelf shaker support system without duplex outlet, each shelf will hold 90 kg. for 3948 and 3962. Must be ordered with duplex outlet (505099 or 505094) |
| 190761* | Three shelf shaker support system includes duplex outlet, each shelf will hold 90 kg for 3948 and 3962 |
| Door Options, factory installed | |
| 190239 | Lexan® Inner Door Kit (five doors) for 3948 and 3962; factory installed |
| 190514 | Door Lock Assembly for 3948 and 3962; factory installed |
| 190597 | Left-Hand Door Swing Kit for 3948; factory installed |
| Stainless Steel Door Glass Covers | |
| 190591 | For 3948 and 3962; factory installed |
| 190892 | For 3948 and 3962** |
| Reinforced Floor Options | |
| 500182*** | Reinforced Floor with Removable Ramp for 3948 and 3962; accommodates a cell roller |
| 190777*** | Same as No. 500182 but factory installed |
| Access Ports, factory installed | |
| 505101 | Port for 3948 and 3907; 38mm I.D.; on center right wall |
| 193004 | 51 mm nominal ID port with cover |
| 190164 | Port for 3962, 61 mm I.D., customer specified location |
| 193005 | 102 mm nominal ID port with cover |
| 193006 | 152 mm nominal ID port with cover |
| Convenience Receptacles, factory installed | |
| 505094 | 230V (includes separate line cord) for 3948 and 3907; located on center rear wall |
| Data Outputs (select one) for 3962, factory installed | |
| 190523 | RS-485 interface (compatible with 1535 Monitor/Alarm only) |
| 190512 | 4-20 milliamp |
| 190543 | 0-5V |
| 190544 | 0-1V |
| Sensaphone Telephone Dialing Systems, interface with standard touch-tone phone | |
| 400047 | For up to four input channels |
| 400134 | For up to eight input channels |

Warranty

We confidently back our environmental chambers with a standard limited two-year parts and labor warranty.

*Rubber isolators replace casters for maximum stability.

**Must be installed by qualified personnel.

***Ramp extends approximately 584mm (23") in front of the chamber.

****Specification sheets for protocol and field validations are available for review upon request.

| Thermo Scientific Cat. No | Description |
|---|---|
| Chart Recorders and Paper | |
| 6 in., 7 Day Chart Recorders for 3948 and 3907; 0°C to 60°C (32°F to 140°F) | |
| 201144 | Single Pen, factory installed |
| 201145 | Single Pen, onsite installation |
| 201146 | Dual Pen, factory installed |
| 201147 | Dual Pen, onsite installation |
| 6 in., 7 Day Chart Recorders for 3962, 0°C to 60°C (32°F to 140°F) | |
| 201156 | Single Pen, 230V, 50/60 Hz |
| 6 in. Chart Paper for Recorders, 50 per box | |
| 180006 | For Single Pen, 0°C to 60°C (32°F to 140°F) |
| 197030 | For Single Pen, -10°C to 70°C (14°F to 158°F) |
| 197075 | For Single or Dual Pen, 0°C to 100°C (32°F to 212°F) |
| RH System Options | |
| 1900391 | Condensate Evaporator for 3948 and 3907; includes separate line cord; 240V |
| 1900139 | Heatless Dryer Kit for 3948 and 3907, for sub-ambient chamber RH, factory installed; The heatless dryer requires an air supply of 6.2 bar (90 PSI) at 17 m3/h (10 CFM) compressor, at minimum. It is recommended that the air supply be greater than the minimum required by the dryer. |
| Deionization Cartridge Assembly for chambers with humidity and no DI water source | |
| 019-168-00 | DI cartridge assembly includes DI cartridge, wall bracket, pressure regulator and gauge and 1/4" flexible tubing |
| 585-036-00 | Replacement DI cartridge |
| Calibration Certificate and Temperature Mapping with Report | |
| 260045 | Temperature Mapping Report (factory installed) |
| 260049 | Calibration Certificate (specify test parameters) |
| cGxP Protocols and on site Validation Services**** | |
| GLP Level Protocols (IQ OQ) | |
| IQQDOCE89003499 | For 3962 |
| IQQDOCE89003710 | For 3948, 3907 |
| GMP Level Protocols (IQ OQ PQ) | |
| IOPQDOCE89003499 | For 3962 |
| IOPQDOCE89003710 | For 3948, 3907 |
| GLP Level Field Validation (IQ OQ) Performed at the customer site (protocol included at no extra charge) | |
| IQQPCKE89003499 | For 3962 |
| IQQPCKE89003710 | For 3948, 3907 |
| GMP Level Field Validation (IQ OQ PQ) Performed at the customer site (protocol included at no extra charge) | |
| IOPQPCKE89003499 | For 3962 |
| IOPQPCKE89003710 | For 3948, 3907 |
| Miscellaneous Accessories | |
| Infrared CO₂ Control Packages, factory installed | |
| 1900227 | For 3948 and 3907 |
| Cell Roll System | |
| 4868 | Three-Tier Cell Roller Base (15 positions), 230V, 50/60 Hz, 757mm W x 706mm H x 620mm F-B (29.8" x 27.8" x 24.4") |
| 190049 | Add-On Tier (5 positions), 29.8"W x 7.1"H x 24.4"F-B (75.7cm x 18.0cm x 62.0cm), customer installed |
| 190777 | Reinforced Floor with Removable Ramp, ramp extends 23.0" (58.4cm), factory installed |
| 500182 | Same as No. 190777 but customer installed |
| 228077 | Rotation Alarm System, includes annunciator jack, factory installed |
| 228078 | Battery Back-Up, provides 24 hours of power if a power failure occurs, factory installed |
| 475560 | 110mm x 285mm Glass Bottles (4 per case) |



thermoscientific.com/chambers

© 2015 Thermo Fisher Scientific Inc. All rights reserved. Sensaphone Company. Lexan is a registered trademark of the General Electric Company. All other trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details.

Worldwide Service and Support

We are committed to keeping your lab equipment working at peak performance levels. Our goal is to help you lower ownership costs, manage labs more effectively, and increase productivity. Contact your Thermo Scientific representative to learn more about our service offerings, including service agreements, preventive maintenance, on-site field repair, depot repair, compliance services and educational services.

North America: USA/Canada +1 866 984 3766 (866-9-THERMO)
www.thermoscientific.com/chambers

Europe:

Austria +43 1 801 40 0,
Belgium +32 53 73 42 41,
France +33 2 2803 2180,
Germany national toll free 08001-536 376,
Germany international +49 6184 90 6940,
Italy +39 02 95059 448,
Netherlands +31 76 579 55 55,
Nordic/Baltic/CIS countries +358 9 329 10200,
Russia +7 812 703 42 15,
Spain/Portugal +34 93 223 09 18,
Switzerland +41 44 454 12 12,
UK/Ireland +44 870 609 9203

Asia:

Australia +61 39757 4300,
China +86 21 6865 4588 or +86 10 8419 3588,
India toll free 1800 22 8374,
India +91 22 6716 2200,
Japan +81 45 453 9220,
New Zealand +64 9 980 6700,
Other Asian countries +852 2885 4613

Countries not listed: +49 6184 90 6940

Thermo
SCIENTIFIC

A Thermo Fisher Scientific Brand